



O2k-Protocols

Mitochondrial Physiology Network 09.12(11): 1 (2014)
Updates: www.bioblast.at/index.php/MiPNet09.12_O2k-Titrations

©2004-2014 OROBOROS®
Version 11: 2014-04-15

Oxygraph-2k manual titrations: SUIT protocols with mitochondrial preparations



OROBOROS INSTRUMENTS Corp, high-resolution respirometry
Schöpfstr 18, A-6020 Innsbruck, Austria
erich.gnaiger@oroboros.at; www.oroboros.at

Oxygraph-2k chamber volume: 2.0 ml

Substrates	Event	Concentration in syringe (solvent)	Storage [°C]	Final conc. in 2 ml	Titration [µl]	Syringe [µl]
Pyruvate	P	2 M (H ₂ O)	fresh	5 mM	5	25
Malate	M	0.4 M (H ₂ O)	-20	0.5 mM	2.5	25
Glutamate	G	2 M (H ₂ O)	-20	10 mM	10	25
Succinate	S	1 M (H ₂ O)	-20	10 mM	20	50
Octanoyl carnitine	Oct	0.1 M (H ₂ O)	-20	0.2 mM	4	10
Ascorbate	As	0.8 M (H ₂ O)	-20	2 mM	5	25
TMPD	Tm	0.2 M (H ₂ O)	-20	0.5 mM	5	25
Cyt. c	c	4 mM (H ₂ O)	-20	10 µM	5	25
ADP+ Mg ²⁺	D	0.5 M (H ₂ O)	-80	1 - 5 mM	4 - 20	25
ATP+ Mg ²⁺	T	0.5 M (H ₂ O)	-80	1 - 5 mM	4 - 20	25
Uncoupler						
CCCP ¹	U	0.1 mM (EtOH)	-20	0.05 µM steps	1 µl steps	10
Inhibitors						
Rotenone	Rot	1.0 or 0.2 mM (EtOH)	-20	0.5 or 0.1 µM	1	10
Malonic acid	Mna	2 M (H ₂ O)	fresh	5 mM	5	25
Antimycin A	Ama	5 mM (EtOH)	-20	2.5 µM	1	10
Myxothiazol	Myx	1 mM (EtOH)	-20	0.5 µM	1	10
Sodium azide	Azd	4 M (H ₂ O)	-20	≥100 mM	≥50	50
KCN	Kcn	1 M (H ₂ O)	fresh	1.0 mM	2	10
Oligomycin	Omy	4 mg/ml (EtOH)	-20	2 µg/ml	1	10
Atractyloside	Atr	50 mM (H ₂ O)	-20	0.75 mM	30	50
Cell perm.						
Digitonin	Dig	10 mg/ml (DMSO)	-20	10 µg·10 ⁻⁶ cells	1µl 10 ⁻⁶	10
Other						
Catalase	Ctl	112,000 U/ml (MiR05)	-20	280 U/ml	5	25

¹ Higher concentrations (1 mM stock) for intact cells in various culture media (e.g. RPMI, DMEM, EGC) and for TIP2k.

Reference: Pesta D, Gnaiger E (2012) High-resolution respirometry. OXPHOS protocols for human cells and permeabilized fibres from small biopsies of human muscle. *Methods Mol Biol* 810: 25-58. »